

Quantitative Measurement of Hexavalent Chromium Treatment to Metal Screws

Introduction

Hexavalent chromium (Cr(VI)) is a significantly hazardous material regulated by OSHA with reference to [29CFR parts 1910 and 1915](#), by RoHS directive, and by [IEC-62321](#) which strictly controls and regulates the use of specific hazardous materials in a wide range of industrial processes as well as electrical and electronic products.

This application note describes the extraction and quantification of residual chromium from the treatment of two samples of screws using a chromogenic reaction and UV-Visible spectroscopy.


The chromium from two steel screws (Figure 1) was extracted via Hydrothermal Extraction. The screws were placed in 25 mL of purified water and heated at 80 degrees Celsius for 30 minutes. These conditions are subject to change depending on the sample type and regulatory compliance.

Keywords

V-630, UV-Visible/NIR, Materials, IEC-62321, 29CFR part 1910, 29CFR part 1915

Experimental

The chromium from two steel screws (Figure 1) was extracted via Hydrothermal Extraction. The screws were placed in 25 mL of purified water and heated at 80 degrees Celsius for 30 minutes. These conditions are subject to change depending on the chromium coatings for a particular screw.

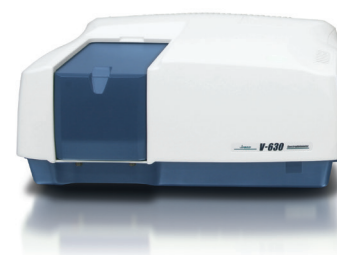

V-630
UV-Visible Spectrophotometer



Figure 1. Sample screws.

After hydrothermal extraction, the screws were removed from the extraction liquid which was cooled to room temperature. The chromogenic reagent (Kyoritsu Chemical-Check Lab) was added to the extraction liquid, stirred for 1 minute, and the left to stand for 5 minutes. Turbidity was observed after 5 minutes. Figure 2 shows the sample pre- and post-filtration. Measurements were obtained on both samples for comparison using the following measurements parameters.

Measurement Conditions			
Bandwidth	1.5 nm	Data Interval	0.2 nm
Scan Speed	400 nm/min	Response	Medium



Figure 2. Screw samples after the chromogenic agent was added to the extraction solutions and pre- (left) filtration and post (right) filtration of the solutions.

Results

The spectra of the samples before and after filtration are illustrated in Figure 3. The spectrum of the sample before filtration shows a higher baseline, which is due to the dispersion of interfering substances in the solution. A three wavelength quantitative analysis was performed to confirm the effects of the interfering substances. Table 1 shows the results of the peak height calculations. Regardless of the pretreatment, almost identical results were obtained.

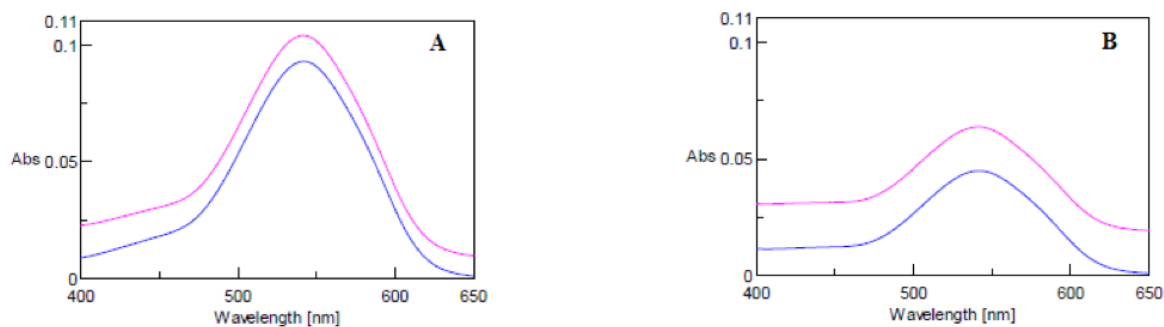


Figure 3. Absorption spectra of samples A and B before (purple) and after (blue) filtration.

Table 1. Peak height calculation results of screw samples A and B.

		Sample A		Sample B	
		Before	After	Before	After
1 Wavelength	Abs at 542 nm	0.10393	0.09288	0.06365	0.04482
3 Wavelength	Peak height: 542 nm Base: 402, 635 nm	0.08782	0.08765	0.03932	0.03915

Three wavelength quantitative analysis was examined on the peak height calculated with peak wavelength at 402 nm and a base wavelength at 635 and 542 nm. The calibration curve obtained by the diphenylcarbazide method in Application Note UV-0004 was applied to this quantitative analysis. Table 2 shows the concentration of Cr (VI) for samples A and B using a two-sided 95% confidence interval and both samples A and B fall within ± 0.005 mg/L of the confidence interval.

Table 2. Quantitative results of Cr(VI) in screw samples.

Sample A		Sample B	
Before	After	Before	After
0.135 \pm 0.006 mg/L	0.135 \pm 0.006 mg/L	0.059 \pm 0.005 mg/L	0.059 \pm 0.005 mg/L